**Lovibond**

# MINIKIT

- Ⓜ **GB Cyanuric Acid**
- Ⓜ **DE Cyanursäure**
- Ⓜ **FR Acide Cyanurique**
- Ⓜ **IT Acido Cianurico**
- Ⓜ **ES Ácido Cianúrico**
- Ⓜ **PT Ácido Cianúrico**
- Ⓜ **DK Cyanursyre**
- Ⓜ **NL Cyanuurzuur**

20 - 200 mg/l

**AF 422**

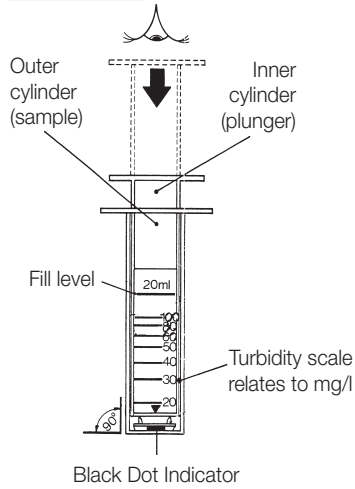
**41 42 20**

**GB**

Assembled unit


**Introduction**

Iso cyanuric acid is used, generally as a constituent of chlorine donors recommended for use in outdoor pools, to stabilize chlorine against loss by the effect of sunlight. It is essential to monitor the cyanuric acid level as too little results in rapid loss of chlorine and too much results in a reduction in the effectiveness of the chlorine as a disinfectant.

**Contents**

- 1 double tube assembly
- 1 stirring rod
- 1 cleaning brush
- 50 CyA-TEST tablets
- 1 operating manual in 8 languages

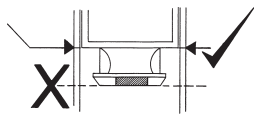**Take reading from shoulder**

## Instruction

1. Separate the double tube assembly.
2. Fill outer cylinder to the 20 ml mark with the water sample.
3. Add a CyA-TEST tablet. Crush the tablet with a clean stirring rod and mix well.
4. Particles which settle at the bottom after the tablet has dissolved can be disregarded. A uniform cloudiness of the water sample indicates the presence of cyanuric acid.
5. Insert the inner tube slowly into the outer cylinder and allow it to fill with the water sample through the two holes at the bottom.
6. Hold the outer cylinder up as far as possible (see illustration) to avoid light losses. The observer looks vertically from above the unit and moves the inner tube gently up and down until the black spot on the base just disappears. Hold the unit so that position of the inner tube is maintained.
7. Read the graduation mark on the outer tube corresponding to the position of the lower edge of the inner tube. This gives the cyanuric acid concentration in the water sample. If the edge is between graduations it is possible to make an estimate of the concentration.

## Notes

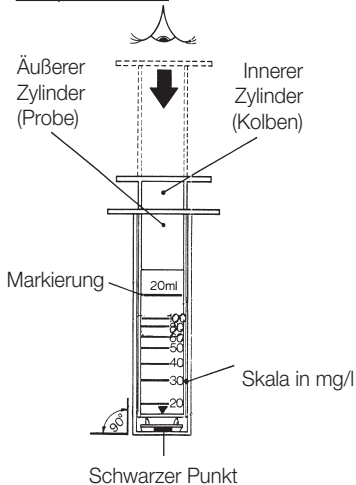
1. After each measurement, thoroughly clean the outer cylinder, flask and stirring rod with the brush included in the kit.
2. If the solution is too cloudy to obtain a reading the concentration is above 100 mg/l cyanuric acid, and the test should be repeated with a diluted water sample. Fill the tube only to the lower (10ml) line. Make up to the top line with cyanuric acid free water (e.g. tap water) and repeat the test as above but multiply the result by 2.

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are Trademarks of the
The Tintometer Ltd.
Tintometer GmbH