LOVIBOND
MINIKIT

GB: Sulphite E: Sulfito
D: Sulfit P: Sulfito
F: Sulfite DK: Sulfit
I : Solfito NL: Sulfiet

20 - 500 mg/l Na₂SO₃

AF 435
41 43 50

GB

Introduction

The Sulphite High Range MINIKIT provides a simple means of checking sulphite levels in water over the range 20 - 500 mg/l sulphite as Na₂SO₃, and it is based on the tablet count method.

Contents

calibrated sample container, 100 ml
Sulfite No.1 tablets, 40 pieces in foil
Sulfite No.2 HR tablets, 100 pieces in foil
operating manual in 8 languages

Instructions

Range 20 - 500 mg/l	Probenvolumen 50 ml
20 - 250 mg/l	100 ml

1. Select the sample size 50 ml or 100 ml appropriate to the range under test. Take a sample of the correct size in the sample container.
2. Add two Sulfite No.1 tablets, stopper the container and shake until the tablets have disintegrated.
3. Add one Sulfite No.2 HR tablet and swirl the container gently until the tablet disintegrates.
Continue adding Sulfite No.2 HR tablets in this manner one at a time until a **BLUE** colour appears.
4. Note the number of Sulfite No.2 HR tablets used and calculate the result from the formula below:

For a 50 ml Sample Size:

Sulphite (as Na₂SO₃) = (No. of Sulfite No.2 HR tablets x 40) - 20

For a 100 ml Sample Size:

Sulphite (as Na₂SO₃) = (No. of Sulfite No.2 HR tablets x 20) - 10

Cleaning

Thoroughly rinse out the sample container after each use.

Example

No. of Sulfite No.2 HR tablets used = 6

Sample Size: 50 ml

(6 x 40) - 20 = 220 mg/l Na₂SO₃

Notes

1. Note the Sulfite No.1 tablets are to condition the sample only. Do not include these tablets in the calculation of sulphite level.
2. For the determination of low levels of sulphite in water use the Sulphite Low Range MINIKIT AF 434. This covers the range 2 - 50 mg/l sulphite as Na₂SO₃.

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