Lovibond® Water Testing

Tintometer® Group

SensoDirect 150


Data Logger/Acquisition Software

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Seite 2–5

GB Instruction Manual

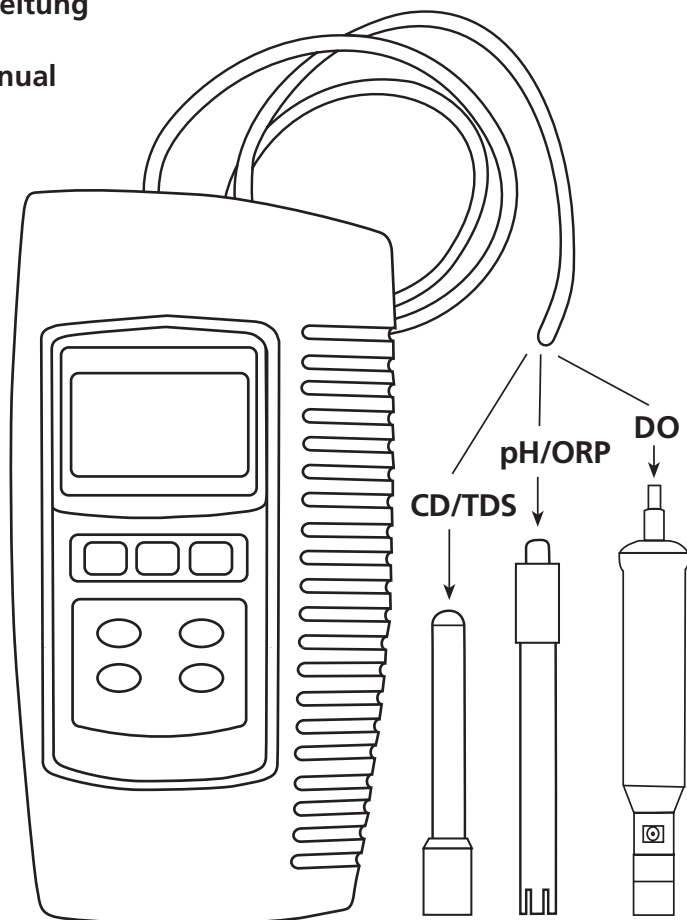Page 6–8

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1.	Allgemeine Beschreibung.....	3
2.	Softwareinstallation	3
3.	SW-E802.xls	3
4.	Meter_Test	5

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
The screenshot shows an Excel spreadsheet with columns A through K and rows 1 through 6. The interface includes several buttons: 'Config / Start' in cell A1, 'stop' in B1, 'Manual Inout' in C1, and 'Clear' in E1. Below these buttons, the spreadsheet has columns for 'Date' (B2), 'Time' (C2), 'Ch1_Value' (D2), 'Ch1_Unit' (E2), 'Ch2_Value' (F2), 'Ch2_Unit' (G2), 'Ch3_Value' (H2), 'Ch3_Unit' (I2), 'Ch4_Value' (J2), and 'Ch4_Unit' (K2). The 'Ch' column is partially visible in column L. The spreadsheet is currently empty, with only the headers and buttons populated.

	A	B	C	D	E	F	G	H	I	J	K	L
1	Config / Start	stop	Manual Inout	Clear								
2		Date	Time	Ch1_Value	Ch1_Unit	Ch2_Value	Ch2_Unit	Ch3_Value	Ch3_Unit	Ch4_Value	Ch4_Unit	Ch
3												
4												
5												
6												

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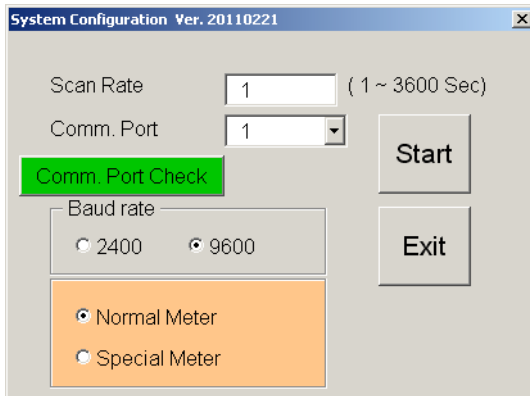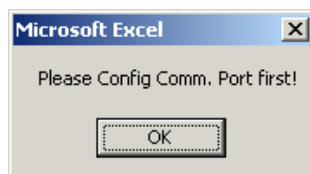
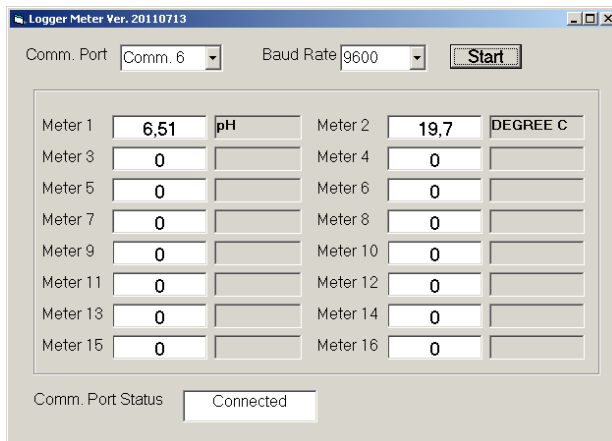
1. General Description

The SW-E802 application software is designed to transfer and log the data generated by the SensoDirect 150 meter in an Excel Spread Sheet. Also an on-line display of the instrument's data on a PC is possible.

An RS232 cable is required to transfer the data. Should the computer not have an RS232 interface, the instrument also can be connected via a USB cable. A virtual COM port can be established using the drivers supplied on the CD.

Part No.: 724500 (RS232 cable)

Part No.: 724510 (USB cable with virtual COM port)

2. Software Installation

If an older version of the software has previously been installed, please delete it before installing the new version. To install the software, insert the CD into the CD ROM disk drive of the computer. After the automatic start-up, please follow the instructions of the installation wizard.


On successful installation, a new programme SW-E802 will appear the "Start/All programs" directory.

3. SW-E802.xls

On execution of the SW-E802.xls, the following Excel-Sheet will appear on the screen.

	A	B	C	D	E	F	G	H	I	J	K
	Confia / Start	Stop Date	Manual Input Time	Ch1_Value	Clear Ch1_Unit	Ch2_Value	Ch2_Unit	Ch3_Value	Ch3_Unit	Ch4_Value	Ch4_Unit
1											
2											
3											
4											
5											
6											

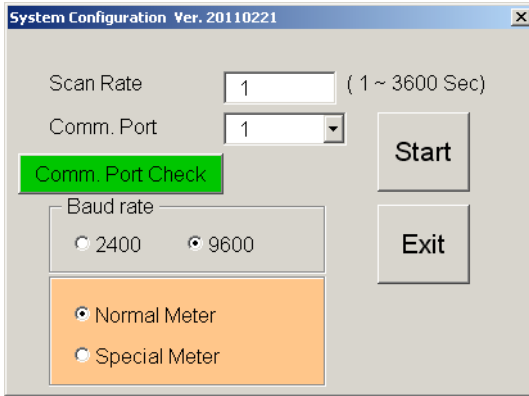Click the 'Config/Start' button to access the below screen where all the required settings can be selected. Confirm the settings with the 'Start' button. The data logging will now begin.


The following settings may be selected:

- Select the scan rate.
- Choose the COM port used to connect the hand held meter. 'Comm. Port Check' enables the user to select the required COM interface.
- Set the baud rate (recommended: 9600).
- Select the meter (Normal meter).
- Select the 'Exit' button to close the setting window without saving the changes.
- Click the 'Start' button to save the settings and start the data logging.

	A	B	C	D	E	F	G	H	I	J	K	
	Config / Start	Stop Date	Manual Input Time	Clear Ch1_Value	Ch1_Unit	Ch2_Value	Ch2_Unit	Ch3_Value	Ch3_Unit	Ch4_Value	Ch4_Unit	Ch
1												
2	1	05.12.2011	08:54:53	5,58	pH	19,7	DEGREE C					
3	2	05.12.2011	08:54:58	5,58	pH	19,7	DEGREE C					
4	3	05.12.2011	08:55:03	5,58	pH	19,7	DEGREE C					
5												
6												
7												

The data will continue to be logged until the 'Stop' button is selected.

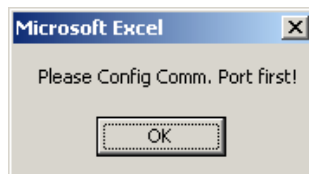The 'Manual Input' button enables the user to store single readings to the Excel-sheet. Transmission is only possible if the settings have been correctly selected. Otherwise the following message appears.


The 'Clear' button can be selected to delete all data logged on the Excel-Sheet. The following warning window will appear. Select 'Ja' to confirm deletion of data or 'Nein' to cancel.

4. Meter_Test

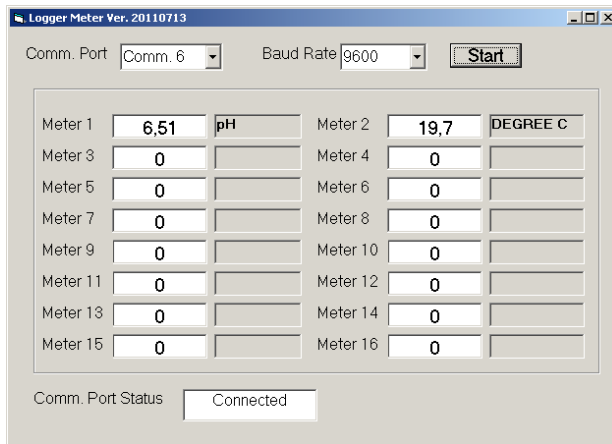On starting Meter_Test, the following operating window is displayed. Once the correct COM port for connection and the baud rate (recommended: 9600) have been selected, the reading of the hand held meter can be displayed by clicking the 'Start' button.

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Technical changes without notice
Printed in Germany 12/11

Lovibond® and Tintometer®
are Trademarks of the
Tintometer Group of Companies

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The Tintometer Limited
Tintometer AG
5212 Hausen AG